



# DeltaVision Ultra


## Switching ON the system

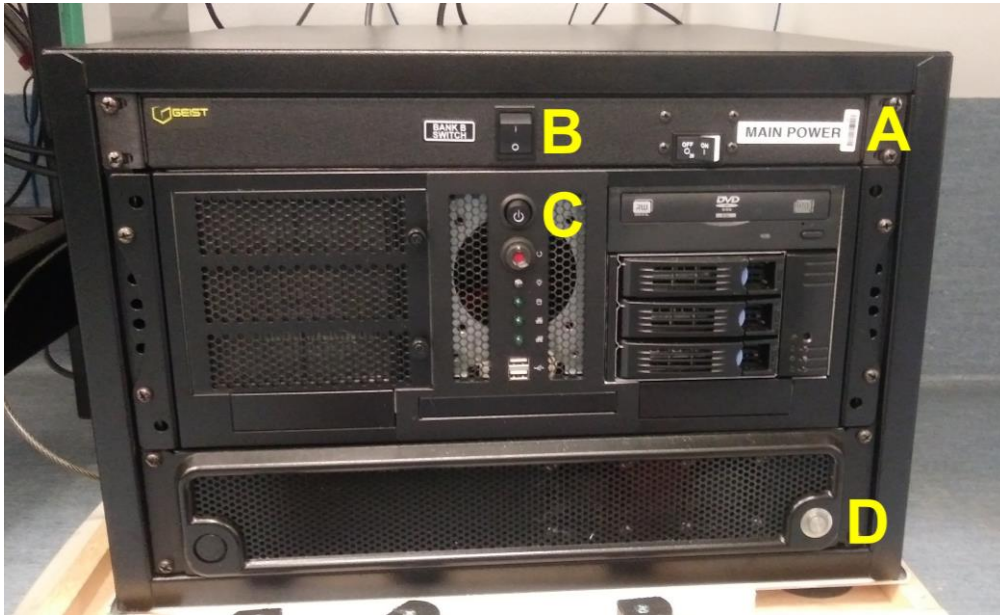

- Switch on the **Main Power** switch (A).
- Switch on the **Bank B** switch (B)
- Switch on the **Workstation Power** switch (C).
- Switch on **M.I.C. (Microscope Interface Chassis)** switch (D). , follow the screen prompts and wait Microscope initialization.
- Log onto the linux CentOS system with your **ALEMBIC Personal Username&Password**.

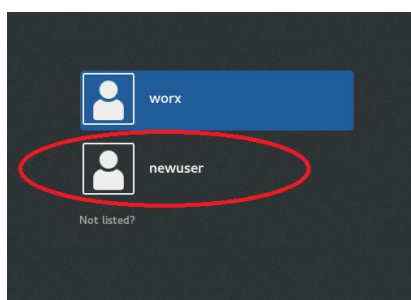

Click the gear (settings) icon near the Sign In button and select **KDE Plasma Workspace**

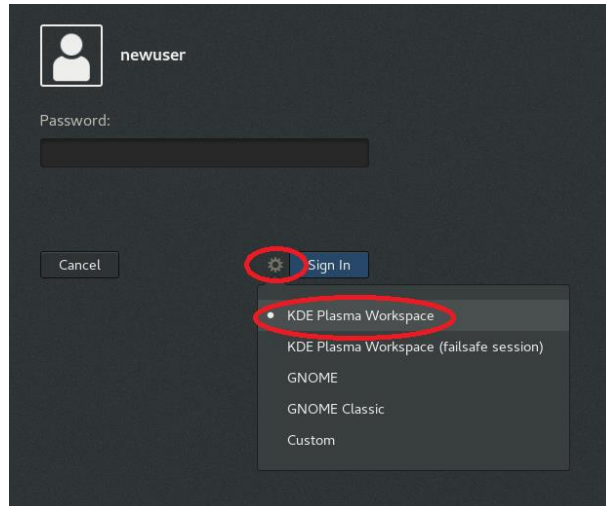

Launch **softWoRx** and **Acquire Ultra** softwares with desktop icons



## Switching OFF the system

If you are the last booked user for the day...

- ✦ Exit from the **Acquire Ultra** and **SotfWorx** softwares
- ✦ Switch off the **M.I.C. (Microscope Interface Chassis)** switch (**D**).
- ✦ Switch off the **Workstation Power** switch (**C**)
- ✦ Switch off the **Bank Bank** switch (**B**)
- ✦ Switch off the **Main Power** switch (**A**).
- ✦ Clean and low the **objectives**.

### ATTENTION:

use clean lens tissue and blot (do not wipe laterally) the lens!




### Tips:

- ✦ All files need to be removed from the DV acquisition computer.  
Store files in the DV Workstation only.

To avoid exhausting the available disk space, remove data as soon as possible.

You can archive your data by writing CD/DVD-ROMs or transfer it via FTP or the Microsoft network.

You can setup a quick connection on Network folder browser

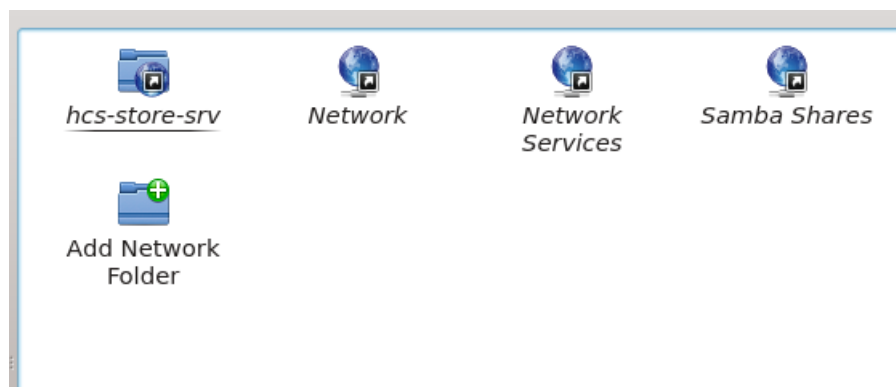

- The **DIBIT FileSharing server**
  - o IP: 172.17.253.30
  - o **srnasric01**
    - and then then select your share space (ex. share01)

Username: hsr\lastname.firstname (*hsr mail's credentials*)

Password: owner

- **ALEMBIC server (hcs-store-srv)**
  - o IP: 172.21.51.90
  - o **hcs-store-srv**

Data on the alembic server are available on any Microsoft, Apple and Unix systems (with ALEMBIC personal Username&Password) and will be automatically deleted after 15 days.