

Mavig RS-G4

Switching ON the system

- Switch on the **Laser Box** (move key clockwise)



N.B.:

Waiting the acoustic signal of system ready before to launch acquisition software

- Switch on the **Prior Proscan III controller** (on the back right of the box)



- Switch on the **PC**
- Log onto the Win 10 system with your **Personal Username and Password**.



Each user should utilize a directory **Data disk**, such as the following:

D:\Data\group name\


- Start **I.D. Software** using the icon on the desktop.



Switching OFF the system

If you are the last booked user for the day...

- Exit from the I.D Software
 - Log-off and close your user session.
- Shut down the computer.

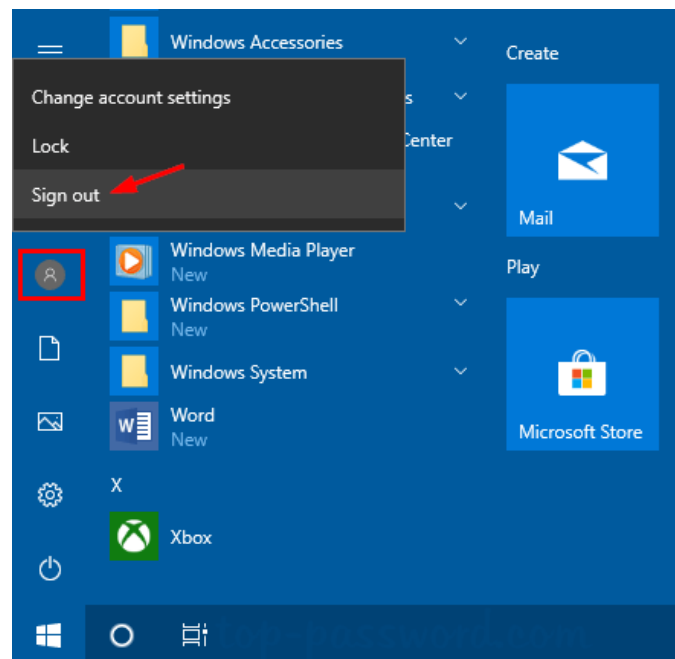

- Remove any objective from the scan head
- Switch off **Prior Proscan III controller** and **Laser Box** (move key counterclockwise)



Tips:

- All images are already saved on the workstation during acquisition. Before finishing your session take care to have time to transfer all your data before the following user

To avoid exhausting the available disk space, remove data as soon as possible. You can archive your data by writing CD/DVD-ROMs or transfer it via FTP or the Microsoft network.

- The **DIBIT FileSharing server**
 - o IP: 172.17.253.30

Type in Menù: **Start -> Run** or in a **folder's address bar**:

\\srnasric01 and then then select your share space (ex. share01)

Username: hsr\lastname.firstname (*hsr mail's credentials*)

Password: owner

- **ALEMBIC server (hcs-store-srv)**
 - o IP: 172.21.51.90

Microsoft Windows Network → hcs-store-srv

or

Type in Menù: **Start -> Run** or in a **folder's address bar**:

\\hcs-store-srv

Data on the alembic server are accessible on any Microsoft, Apple and Unix systems (with ALEMBIC personal Username&Password) and will be automatically deleted after 15 days.

.